

CALCIUM INDICATORS SELECTION GUIDE

Indicator Dye	MW	Ex	Em	Laser	Filter	Kd	FCa/FFree*
Cal-500™	1019.92	388	482	405	DAPI	303 nM	~100 fold
Calbryte™ 520	1090.9	493	515	488	FITC	1.2 μM	~300 fold
Cal-520®	1102.95	493	515	488	FITC	320 nM	~100 fold
Calbryte™-520L	794.88	493	515	488	FITC	91 μM	~300 fold
Calbryte™-520XL	1129.36	493	515	488	FITC	300 μM	~300 fold
Cal-520FF™	1138.92	493	515	488	FITC	9.8 μM	~100 fold
Cal-520N™	1147.96	493	515	488	FITC	90 μM	~100 fold
OG488 BAPTA -1	1258.07	493	522	488	FITC	170 nM	~14 fold
Fluo-8®	1046.93	495	516	488	FITC	389 nM	~200 fold
Fluo-8H™	1074.98	495	516	488	FITC	232 nM	~200 fold
Fluo-8L™	1078.95	495	516	488	FITC	1.9 μM	~200 fold
Fluo-8FF™	1082.91	495	516	488	FITC	10 μM	~200 fold
Fluo-4	1096.95	495	528	488	FITC	345 nM	~100 fold
Cal Green™ 1	1290.96	498	517	488	FITC	190 nM	~14 fold
Fluo-3	1129.85	506	515	488	FITC	390 nM	~100 fold
Rhod-4™	1015.96	523	551	561	TRITC	451 nM	~250 fold
Rhod-2	1123.96	553	577	561	TRITC	570 nM	~100 fold
Rhod-FF	1145.9	553	577	561	TRITC	19 μM	~100 fold
Rhod-5N	1154.92	557	580	561	TRITC	300 μM	~100 fold
Cal-590™	1266.81	574	588	561	Cy3/TRITC	561 nM	~100 fold
Calbryte™ 590	1218.77	581	593	561	Cy3/TRITC	1.4 μM	~300 fold
Calbryte™ 630	1234.84	607	624	561	Texas Red	1.2 μM	~300 fold
Cal-630™	1282.89	609	626	561	Texas Red	792 nM	~100 fold
Cal-670™	1587.99	667	680	640	Cy5	853 nM	~100 fold
Cal-770™	1688.58	758	783	750	Cy7	850 nM	~100 fold

* FCa/FFree = Increase in fluorescence intensity of calcium bound indicator relative to its calcium free indicator form.

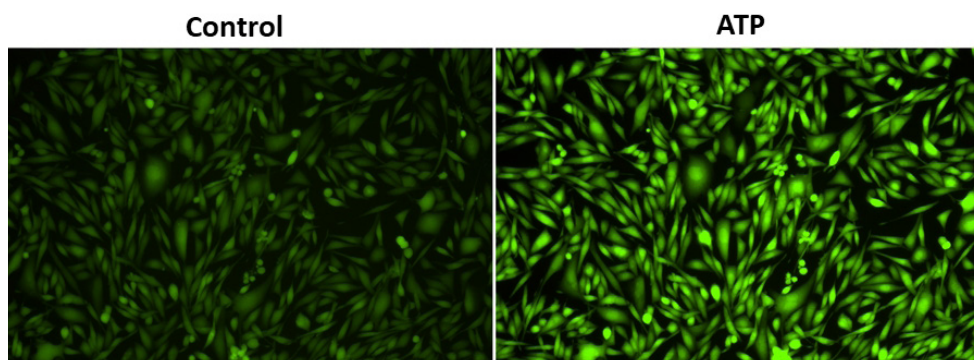


Figure 1. Response of endogenous P2Y receptor to ATP in CHO-K cells. CHO-K cells were seeded overnight at 40,000 cells per 100 μL per well in a 96-well black wall/clear bottom costar plate. 100 μL of 4 μM Cal 520™ AM in HHBS with 1 mM probenecid were added into the wells, and the cells were incubated at 37 °C for 1 hour. The dye loading mediums were replaced with 100 μL HHBS and 1 mM probenecid, then imaged with a fluorescence microscope (Olympus IX71) using FITC channel before and after adding 50 μL of 300 μM ATP.

For Research Use Only. Not for use in diagnostic or therapeutic procedures. © 2024 AAT Bioquest, Inc. All rights reserved. All trademarks and registered trademarks mentioned herein are the property of their respective owners. All other trademarks and registered trademarks are the property of AAT Bioquest.

